


Stable Fluorinating Reagent with Ease of Handling FLUOLEAD™

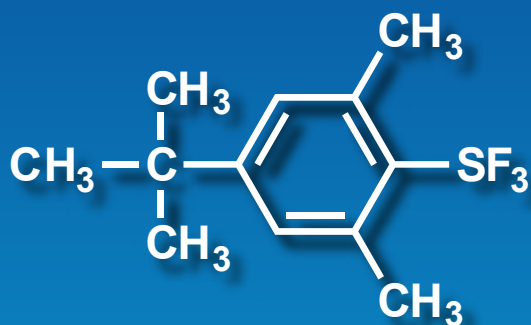

FLUOLEAD™

1g / 5g / 25g

[B3664]

Advantages

- Thermally Stable
- Easy to Handle
- Widely Applicable to Various Substrates

4-*tert*-Butyl-2,6-dimethylphenylsulfur Trifluoride (FLUOLEAD™) [B3664], which was first reported by Umemoto, is a novel nucleophilic fluorinating reagent. Differing from other existing fluorinating reagents, such as DAST, B3664 is a crystalline solid with high thermal stability, which makes it easier to handle. B3664 fluorinates a variety of substrates to afford the corresponding fluorinated compounds in good yields.


Application

Nucleophilic Fluorinating Reagent

- | | | | | | |
|----------|---|-------------------|--|---|---|
| • R-OH | → | R-F | • R-CHO | → | R-CF ₂ H |
| • R-COOH | → | R-CF ₃ | • R ¹ -C(=S)-R ² | → | R ¹ -CF ₂ -R ² |

Example 1


Fluorination of Alcohols¹⁾

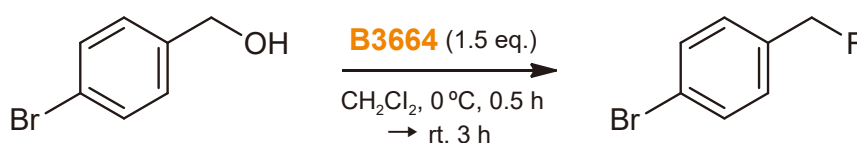

Stable Fluorinating Reagent with Ease of Handling: FLUOLEAD™


Example 2

Fluorination of Aldehydes¹⁾

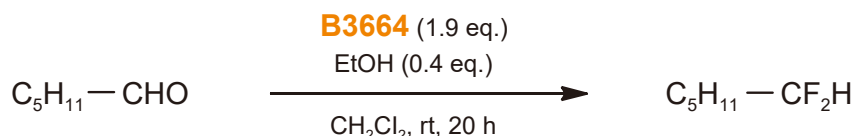

Example 3

Direct Conversion of Carboxyl Group to Trifluoromethyl Group¹⁾

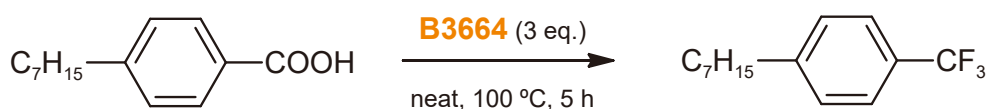

Application Example

Fluorination of *p*-Bromobenzyl Alcohol¹⁾

Into a solution of 7.51 g (30 mmol) of **B3664** in dry dichloromethane (20 mL) cooled on an ice bath, is slowly added a solution of 3.74 g (20 mmol) of *p*-bromobenzyl alcohol in 15 mL of dry dichloromethane. After 0.5 hour, the ice bath is removed and the reaction mixture is stirred at room temperature for 3 hours. The reaction mixture is poured into cooled 100 mL of 5% NaOH aqueous solution and the mixture is stirred for 1 hour at room temperature (note: the byproduct, ArS(O)F can be removed by this alkaline treatment). The product is extracted with dichloromethane and washed with water. The removal of the solvent gives the crude product, which is further purified by chromatography to give 3.0 g of pure *p*-bromobenzyl fluoride (yield: 80%).

For further information please refer to our website at www.TCIchemicals.com. ▶▶▶

TCI fluorination



Ordering and Customer Service

TCI AMERICA

Tel : 800-423-8616 / 503-283-1681
Fax : 888-520-1075 / 503-283-1987
E-mail : Sales-US@TCIchemicals.com

TCI EUROPE N.V.

Tel : +32 (0)3 735 07 00
Fax : +32 (0)3 735 07 01
E-mail : Sales-EU@TCIchemicals.com

TCI Deutschland GmbH

Tel : +49 (0)6196 64053-00
Fax : +49 (0)6196 64053-01
E-mail : Sales-DE@TCIchemicals.com

Tokyo Chemical Industry UK Ltd.

Tel : +44 (0)1865 78 45 60
E-mail : Sales-UK@TCIchemicals.com

梯希爱(上海)化成工业发展有限公司

Tel : 800-988-0390 / 021-67121386
Fax : 021-6712-1385
E-mail : Sales-CN@TCIchemicals.com

Tokyo Chemical Industry (India) Pvt. Ltd.

Tel : 1800 425 7889 / 044-2262 0909
E-mail : Sales-IN@TCIchemicals.com

TOKYO CHEMICAL INDUSTRY CO., LTD.

Tel : +81 (0)3-5640-8878
E-mail : globalbusiness@TCIchemicals.com

* Chemicals itemized in this brochure are for research and testing use only. Please avoid use other than by chemically knowledgeable professionals. * Information such as listed products and its specifications and so on are subject to change without prior notice. * The contents may not be reproduced or duplicated in whole or in part without permission of Tokyo Chemical Industry Co., Ltd.